

APL 416

Applicator for contact plates

BIOMAXIMA presents contact plates applicator for standardized sampling and surface cleanliness testing.

The applicator was prepared for microbiological laboratories, that examine the microbiological purity as required by PN-EN ISO 18593 and pharmacopoeia.

Standardized (500g) pressure and time of sampling (10 seconds) to minimize human error occurring during the execution of test. Credible and repeatable results can be obtained with the applicator, regardless of the type of examination.



1. Applicator off



2. Insert one closed BIOMAXIMA contact plate, then remove the lid



3. Push the applicator to the surface.
LED indicator turns **RED**



4. After 10 seconds LED indicator turns **GREEN**.
Lift the applicator, put the lid back on.